

# Allegro™ Connect Bulk Fill System

# Contents

Introduction	3
The Allegro Connect Bulk Fill System	3
The solution	4
Minimizing non-recoverable volumes of high-value products	13
Scale the equipment to meet your process needs	14
A total bulk fill solution	16
Technical specifications	18
Ordering information	23



# Introduction

Executing any process without the right degree of monitoring, control or reporting, runs the risk of process deviations. This could potentially lead to the adulteration of valuable drug products and intermediates, whilst wasting valuable operator time. Allegro Connect Systems provide robust, accurate and automated platforms that integrate with your manufacturing processes on every level, keeping unit operations within critical parameters to ensure that you spend less time collating data and more time optimizing your process.

Our range of Allegro Connect Systems share a compact form factor, designed with operators in mind, that are configurable to meet your process requirements without the time-consuming pain of modification.

## The Allegro Connect Bulk Fill System

The Allegro Connect Bulk Fill System is an automated filtration and filling system designed to accurately and aseptically aliquot bulk product into multiple biocontainer bags or bottles. Minimizing non-recoverable volume within the flow path is central to the design of the system. Automated techniques and novel design features have been incorporated to ensure that bulk fill step yields are consistent and optimal across a variety of bulk filling scenarios.


Configurable to your needs, the system supports a wide range of liquid filter capsules and various filter combinations. The system can fill up to 320 containers with bulk drug substance, process buffer or cell culture media using preconfigured standard accessories. Designed to minimize process risk, the system dispenses product in a sterile single-use flow path, maintaining a closed process at all times. It automatically vents and wets liquid filters, along with performing filter integrity and manifold leak testing. Process risk is further minimized by a comprehensive validation package characterizing extractables as well as particulates, in accordance with current industry expectations.

Uncontrolled risk and preventable product loss at any stage in the manufacturing process is unwelcome. This is accentuated at this critical stage in the process and highlights the need for an efficient, robust and versatile solution to the bulk filling process.

## The problem

With the cost of manufacturing under constant scrutiny, processes and technologies are consistently being reviewed to increase process productivity and the volume of drug substance produced per year (throughput). Current trends show an increasing number of multi-product facilities being built and/or being retrofitted together with significant investments in automated bioprocessing equipment.

Which of these terms best describes the process at your site for drug substance filling?

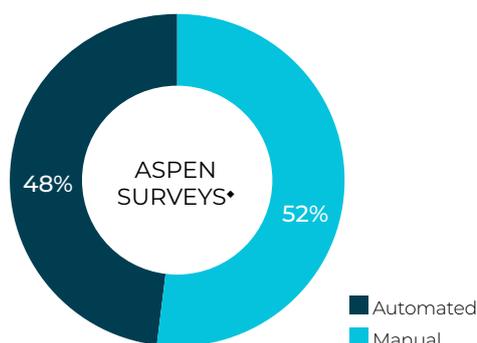

The majority of bulk filling methods currently being used are manual and these potentially limit the production capacity. It is not uncommon to hear of these processes taking up to 30 hours, requiring up to four personnel at any one time. Current automated solutions, however, go some way to address this issue but have single-mindedly focused on throughput of the facility at the expense of step yield and flexibility. A persistent challenge for bulk filling operations is process variability. This is not well addressed by the current generation of automated systems and is one of the main reasons for the relatively low adoption rates of automation in this operation when compared to other bioprocessing equipment.

# The solution

## Minimize non-recoverable volume

Compact flow path

Recovery of product from flow paths before changeout, the filter system and final filter core

Re-dispensing of recovered product into bulk container

## Minimize process risk

Single-use flow path (no change-over) with sterile claim

Processing efficiency gain – minimal human interaction

All data in one place

*In-situ* filter testing and leak testing

## Maximize configurability

Biocontainer bags and bottles (1 L to 20 L)

Up to 320 containers

Choice of manifolds



## The solution is a configurable platform

The Allegro Connect Bulk Fill System is a highly configurable platform that can fill drug substance in both biocontainer bags and bottles. Here are some of the key features the system has to offer:

- Automatically fills up to 320 containers – biocontainer bags or bottles, from 1 L to 20 L nominal volumes.
- Can support single, serial, or parallel liquid filtration arrangements using up to 2 × 508 mm (20 in.) filter capsules with automated venting, priming and aseptic sampling from the flow path via a sampling port upstream of the final liquid filter.
- Non-recoverable product volume within the system is <2.5% of the feed volume.
- If homogeneity of the filtered product is a concern, the system can control an optional mixer, connected aseptically in the flow path after the final filter but before the dispensing station, where a secondary pump is used to dispense the filtered product into the bulk containers.
- It can achieve a filling accuracy of within  $\pm 2\%$ , using a single-floor weighing platform per workstation as opposed to one balance per container. This avoids excessive (re-)calibration activities when a multiple scales system is used for the filling process.

## Dispensing accuracies for containers

Container nominal volume	Biocontainer bags	Bottles
1 L	± 2%	± 2%
2 L	± 2%	± 1%
5 L	± 1%	± 0.5%
10 L	± 0.5%	± 0.3%
20 L	± 0.5%	± 0.3%

- It has an integrated Palltronic® Flowstar IV Integrity Test Instrument to perform pre-use, post-sterilization integrity testing (PUPSIT) and post-use integrity testing of liquid filter capsules. Leak testing of the single-use (SU) flow paths is also carried out, ensuring the main manifold is defect-free.
- It includes a label printer, allowing *in-situ* tagging for all product and sample containers, enabling robust tracing with configurable labels automatically generated.
- It generates electronic batch records along with a summary compiling data from:
  - the single-use sensors within the flow path.
  - single-use manifolds used for the batch.

## Batch record summary for single-use manifold

Component	Part number	Lot number	Serial number	Expiry date
Filter	NP7UECVPIG	IF5156	00012	10-May-2023
Manifold	9496-1356Z	104226303	—	24-Mar-2023
Manifold	9496-1362T	104226077	—	08-May-2023
Manifold	9496-1402Z	104270531	—	10-May-2023
Manifold	7496-1579U	104250435	—	15-Apr-2023
Manifold	7496-1579U	104250435	—	15-Apr-2023
Manifold	9496-1519E	104250572	—	14-Apr-2023

- integrity and manifold leak testing results.

## Integrity test results


Time	Executed test	Message	Unit	Result	Measured value
12-Aug-21 13:17:19	Forward Flow Test – F2 (to Flush)	—	MBF01	Pass	6.5 mL/min
12-Aug-21 15:03:25	Forward Flow Test – F2 (to Flush)	—	MBF01	Pass	5.9 mL/min
12-Aug-21 15:27:05	Leak Test – Up to Primary	—	MBF01	Pass	0.2 mL/min

- batch, lot information and dispensed weight per container, as shown in sample label below.

<b>Campaign ID:</b>	Batch Management	
<b>Description:</b>	Liquid Medicine on	
<b>Fill Date &amp; Time:</b>	DD.MM.YY-09:36	
<b>Container #:</b>	023	
<b>Batch ID:</b>	106	
<b>Storage Temp:</b>	27°C	
<b>Container Tag:</b>	A1 B1 C4	
<b>Lot ID:</b>	102554687	
<b>Fluid Weight:</b>	XXX.XXX kg	
<b>Material #:</b>	Insulin	

- other critical process parameters (e.g., differential pressure across filters).

### Example of the critical process parameter data available

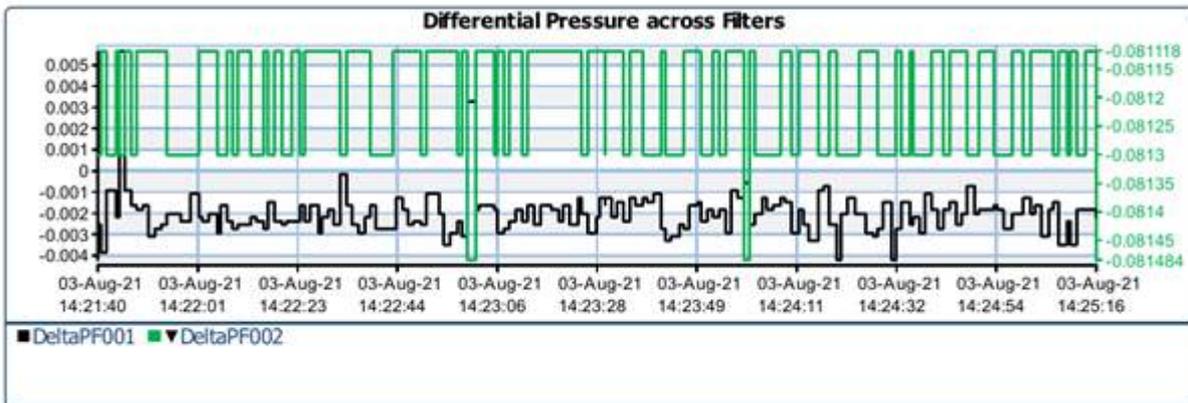

- It controls the container filling flow rate between 0.15 and 10 L/min ensuring (1) the liquid filter capsules are primed correctly, (2) shear damage of the product is avoided, and, if applicable, (3) the processing time for filling the containers is minimized.
- A complete single-use flow path can be installed or uninstalled by a single operator in less than 60 minutes.
- Manual and automated control modes of operation are possible.

# High-performance SU filling

**Processing efficiency gains**  
Up to 80%<sup>1</sup> time saving

80%

**High dispensing accuracies**  
Within  $\pm 2\%$ <sup>2</sup>

$\pm 2\%$

**Minimized product loss**  
< 2.5% of feed volume

< 2.5%

**Compact footprint**  
1.3 m x 2.6 m




<sup>1</sup> Value based on customer feedback. <sup>2</sup> For data refer to Pall Application Note reference USD3669.



## Configure the equipment to meet your process needs


The following images show the flexibility of the system to achieve various configuration options, tailored to meet your individual process needs. With a range of manifolds and hardware configurations available, end users are able to fill up to 320 biocontainer bags and/or bottles per batch.

### Scalable platform for biocontainer bags showing all-in-one and tower systems

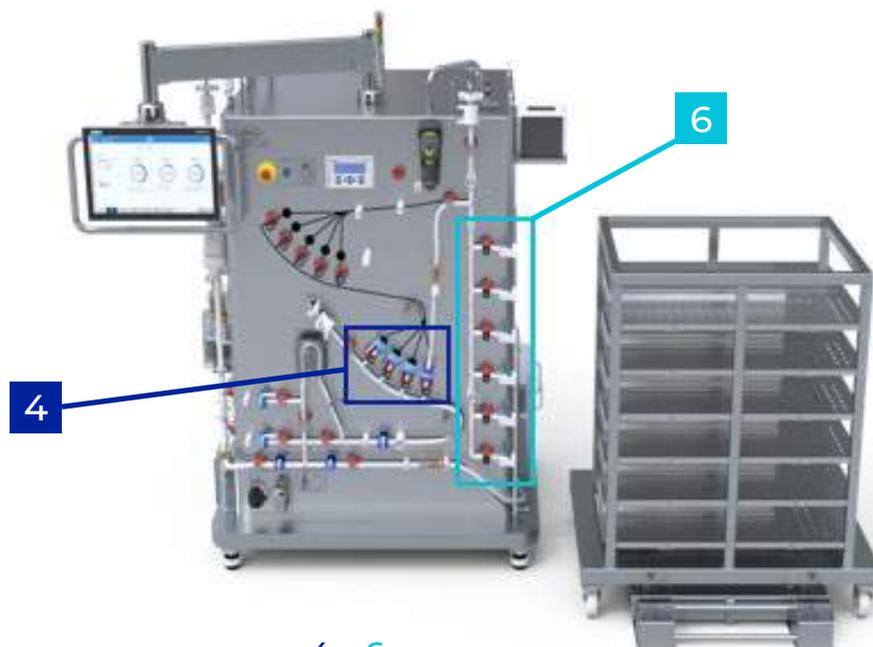


4 x 6

Filling 24 biocontainer bags

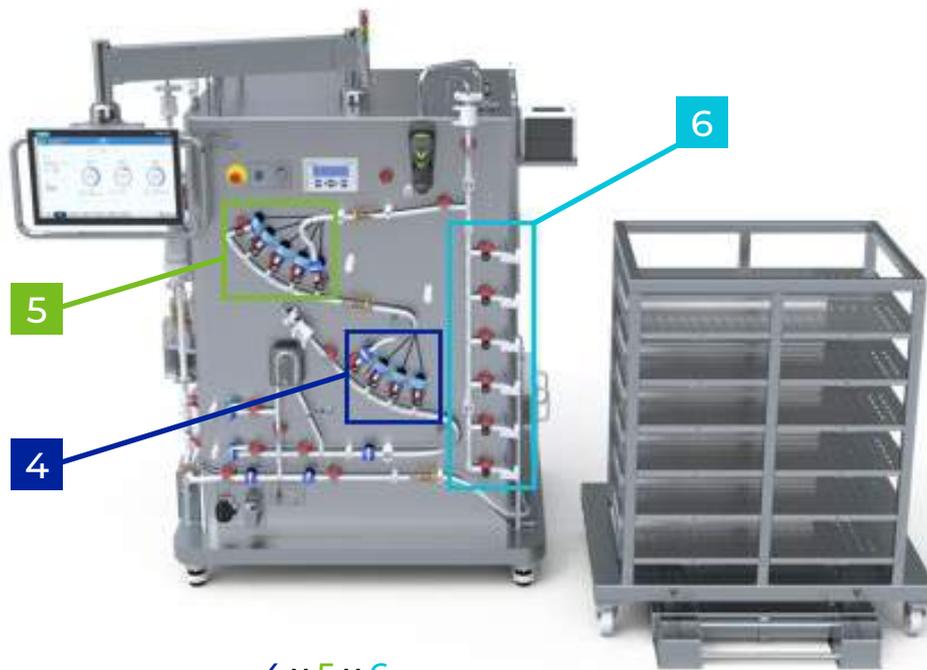



**4 x 5 x 6**  
Filling 120 biocontainer bags


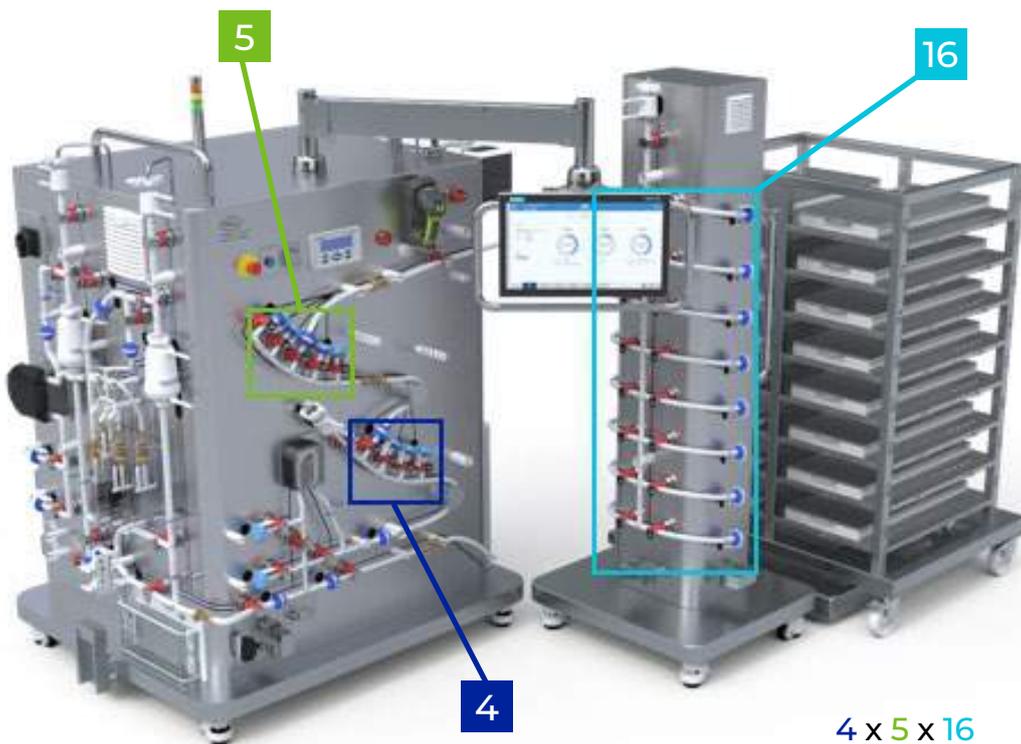


**4 x 5 x 16**  
Filling 320 biocontainer bags\*

\* Additional weighing scale and 16-biocontainer bag workstation required for 10 L and 20 L biocontainer bags

# Scalable platform for bottles showing all-in-one and tower systems

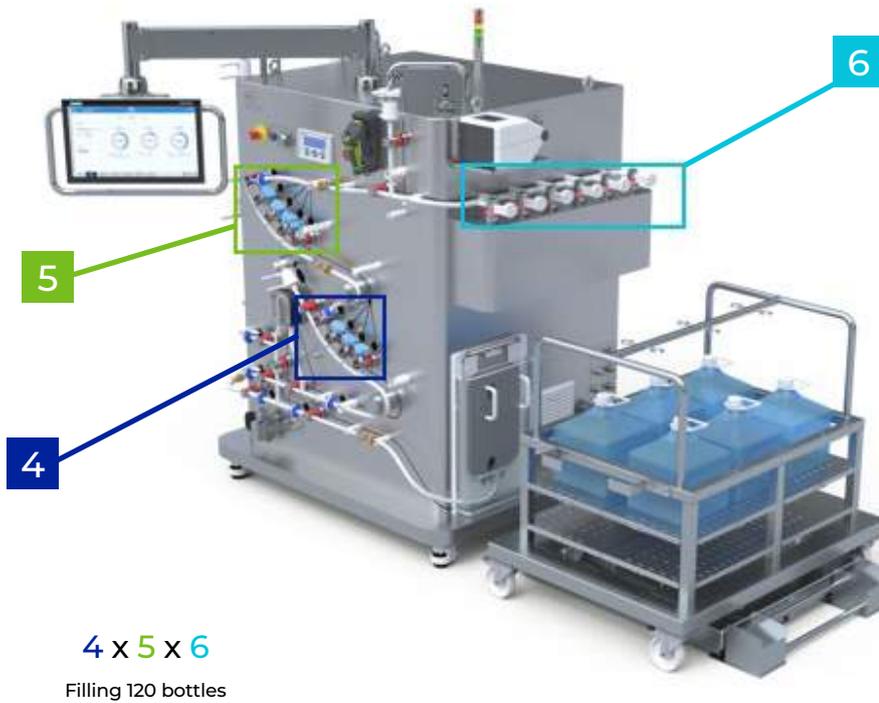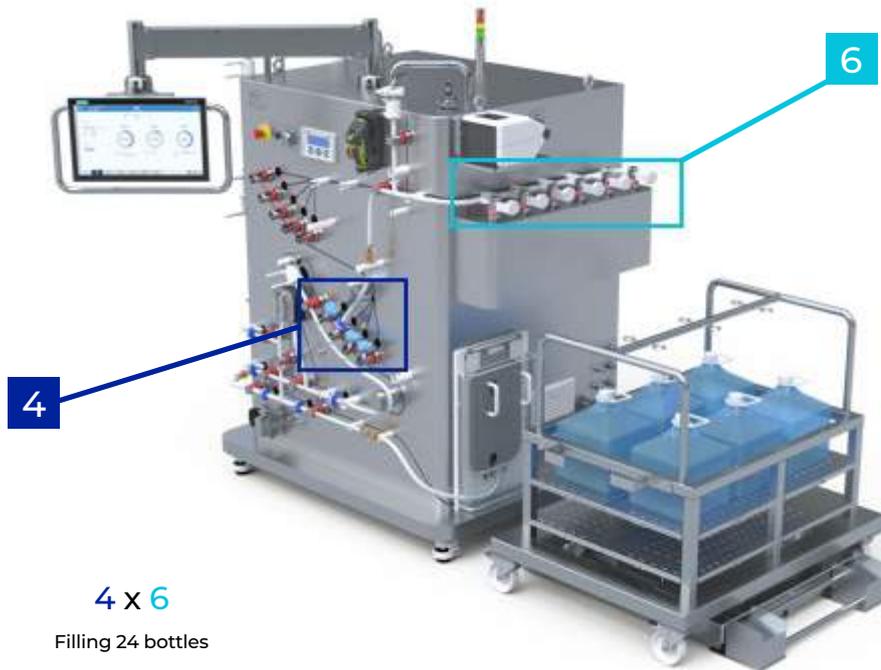

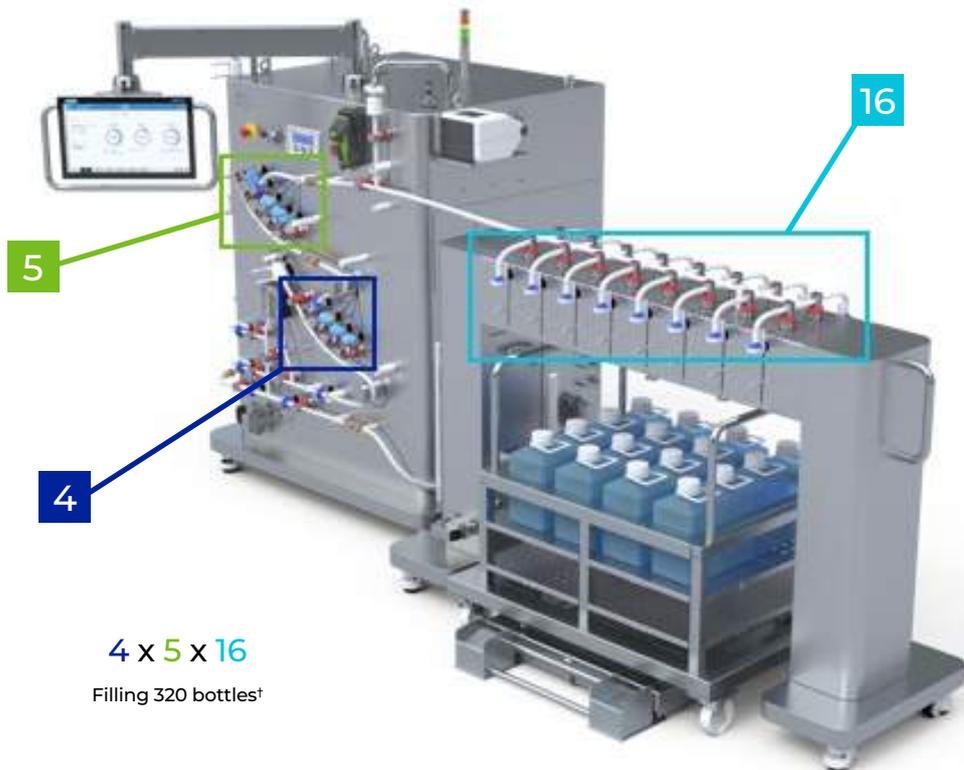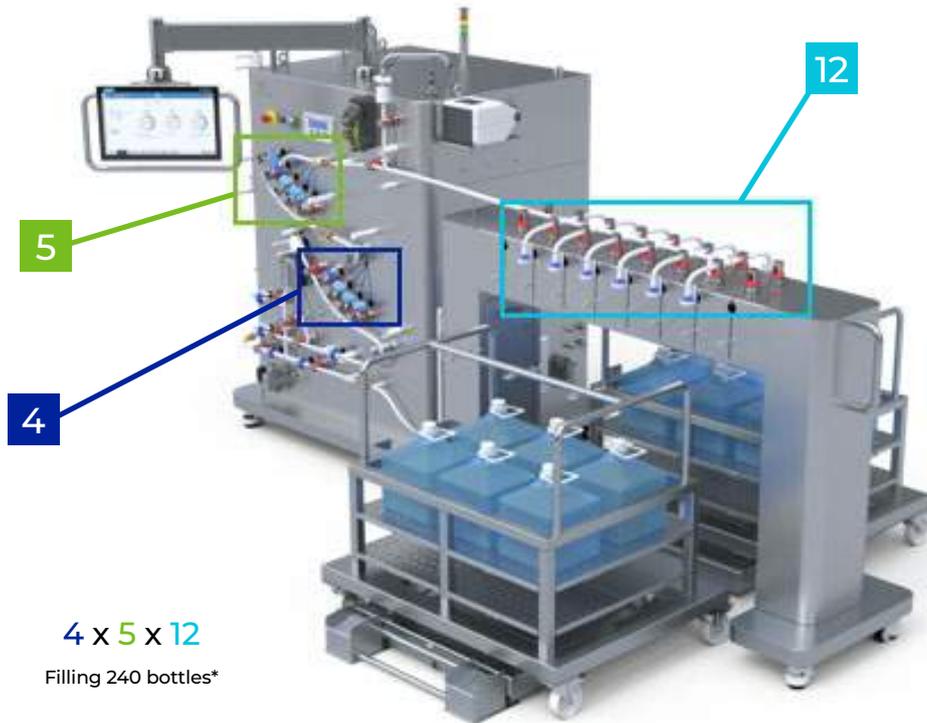

\* Applicable for 10 L and 20 L bottles with 12-bottles workstation (2 trolley + 2 scale).

† Applicable for 1 L to 5 L bottles with 16-bottle workstation (1 trolley + 1 scale).

# Automated manifold leak test and filter integrity test

Automatic leak testing and filter integrity testing completed by the integrated Palltronic Flowstar IV integrity test instrument, with Human Machine Interface (HMI) visualization of the test progress running from pressurization, stabilization, measurement, venting and results.



## Configurable filtration options



No filtration



Single filtration



Serial filtration




Parallel filtration

# Minimizing non-recoverable volumes of high-value products

## A solution you can depend on


### Accuracy

Filling accuracy over a range of container sizes

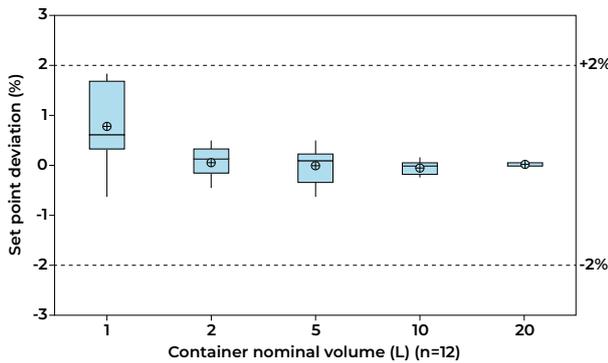

### Consistency

Filling accuracy over varied drug substance viscosity

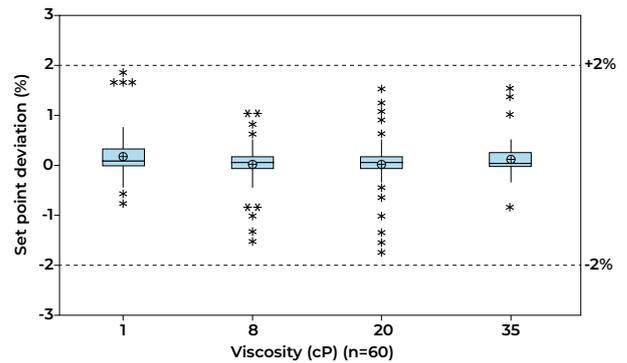


Biocontainer bags and bottles filled to 60% and 80% nominal volume, respectively. The graphs show the accuracy is within  $\pm 2\%$  of the target fill weight for all container sizes and various types of viscous feed stream.

## A solution for all bulk filling scenarios


### Reliability

Non-recoverable product weight by hardware and filtration configuration

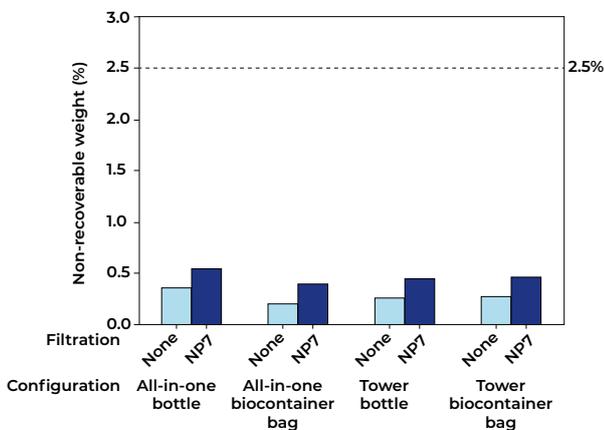

### Range

Non-recoverable product weight by drug substance viscosity

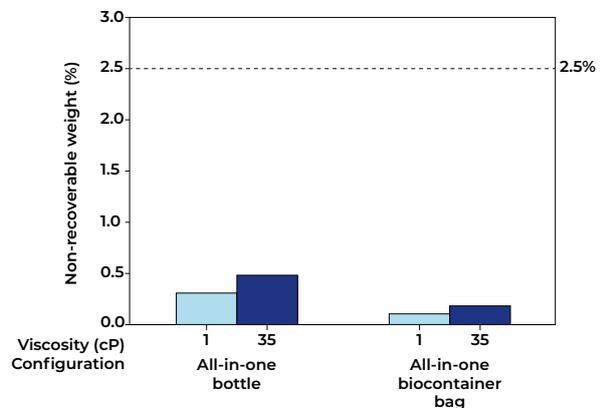

Non-recoverable batch weight shown for a 200 L batch, filling 120 containers with each system set-up and feed type. A no-filtration configuration was used unless otherwise specified. The non-recoverable volume stayed  $<2.5\%$  with different system configuration and with the best-case and worst-case viscosity feed stream.

# Scale the equipment to meet your process needs

## Maximizing productivity

The Allegro Connect Bulk Fill System utilizes single-use technology (SUT) to achieve faster turnaround times between product batches, eliminating the need for clean-in-place (CIP) and sterilization-in-place (SIP) operations and associated cleaning validation. This reduces maintenance cost and system downtime, helping with improved facility throughput.

The entire flow path is designed for easy installation and removal, with clearly marked connections and a shadow board to visibly guide the user and reduce the installation error. Minimal hold-up-volume and optimized flow path design, ensures minimized non-recoverable volumes.



Shadow board concept on the Allegro Connect Bulk Fill System

## Increased assurance

The Allegro Connect Bulk Fill System is supported by a comprehensive validation package including:

- Sterile claim
- Extractables per BPOG
- Particulates requirement per USP <788> aligned with industry expectations (for data refer to regulatory dossier)

Consistency across different batches is achieved through automation and a recipe-driven approach. All batch data is recorded electronically to eliminate any transcription errors.



Allegro Connect Bulk Fill System showing single-use manifolds with Kleenpak® Presto Sterile Connectors and Kleenpak Sterile Disconnectors

## Designed for ease of use

The Allegro Connect Bulk Fill System has undergone extensive usability trials, to ensure the system is simple and intuitive to use.



Guided Instructions for Installation (IFI), where the operator is instructed by the recipe to perform the distribution manifold connection onto the correct valve.



The bulk containers can be assigned to respective valves with the help of the handheld scanner for seamless labeling post filling.



The product container labels are printed with the valve numbers shown in bold font, significantly reducing risk of human error.

# A total bulk fill solution

The Allegro Connect Bulk Fill System can also be connected with some of our other single-use systems to provide a total bulk fill solution for your process, such as the LevMixer® System or Allegro Mixers as a feed tank and intermediate tank, and the Allegro buffer totes for buffer or water for injection (WFI) storage.



## System options

The Allegro Connect Bulk Fill System is available with three types of automation options:

- PLC (Rockwell\* or Siemens\*) and HMI for local stand-alone control
- Remote I/O (no PLC) for integration into a DCS or SCADA system
- Remote I/O (no PLC) controlled by centralized PLC system

The automation architecture is based on either Siemens S7 PLC or Rockwell CompactLogix\* PLC, an industrial PC, and a 22 in thin client which is applied across our range of Allegro Connect bioprocessing systems enabling a truly modular 'plug-and-play' concept, with the ability to control single or multiple unit operations from one centralized cabinet.

The Allegro Connect Bulk Fill System is compatible with all Pall Corporation single-use mixing technology and can also be integrated with most mixing equipment from other suppliers.



## Quality standards

The Allegro Connect Bulk Fill System provides a detailed validation package for all configurations according to ASTM 2500 Standards (a standard guide for specification, design, and verification of pharmaceutical and biopharmaceutical manufacturing systems and equipment).

The regulatory dossier includes:

- Regulatory compliance ROHS I to ROHS III directives
- Raw material compliance data (USP Standards)
- Packaging and packaging waste directive 94/62/EC
- System designed in accordance with the American Society of Civil Engineers (ASCE), ASCE 7, Minimum Design Loads and Associated Criteria for Buildings and Other Structures.

The Pall automation platform enables compliance with 21 CFR Part 11 and follows the GAMP V life cycle for software development.

# Technical specifications

## Process specifications

Equipment	Specification
Functionality	Filtration and/or filling
Filtration configurations	No filter; single; serial or parallel/redundant
Minimum liquid filter capsule size	~ 25 mm (1 in.)
Maximum liquid filter capsule size	~ 508 mm (20 in.)
Number of pumps	2
Pump 1 flow rate range	0.15 – 10 L/min
Pump 2 flow rate range	0.12 – 3.5 L/min
Number of inlets	2 (1 × WFI/buffer and 1 × product)
Number of outlets	Filling: maximum of 16 active outlets, Waste: 2 process waste outlets
Container types to be filled	Biocontainer bags and/or bottles
Tube dimension	Combination of ¼ in. (bottle cap manifold, 1 L and 2 L biocontainer bags), ½ in. and ¾ in. internal diameter (ID)
Flow path operating pressure	0 – 1.5 bar (0 – 22 psi)
Integrity test pressure <sup>1</sup>	Up to 5.7 bar (83 psi) for maximum exposure ≤10 hours
Flow path operating temperature range	4 – 40 °C
Pressure sensor	Up to 4 (1 after each pump with integrated pressure switch; 1 before and after each liquid filter)
External connections	9 (6 × weighing platforms, 2 × mixers, 1 × distribution tower)

<sup>1</sup> Acceptable pressure rating of the air filter

## System specifications

System	Specification
Main system dimensions <sup>2</sup> (W × D × H)	142 × 122 × 199 cm (56 × 48 × 78 in.)
Main system mass (empty)	750 kg (1654 lb)
12-biocontainer bag workstation dimensions (W × D × H)	120 × 90 × 128 cm (47 × 35 × 50 in.)
12-biocontainer bag workstation weight (empty)	165 kg (363 lb)
16-biocontainer bag workstation dimensions (W × D × H)	120 × 90 × 160 cm (47 × 35 × 63 in.)
16-biocontainer bag workstation weight (empty)	220 kg (485 lb)
Bottle workstation dimensions (W × D × H)	120 × 90 × 103 cm (47 × 35 × 40 in.)
Bottle workstation weight (empty)	100 kg (221 lb)
Biocontainer bag distribution tower dimensions (W × D × H)	80 × 70 × 200 cm (31 × 27 × 79 in.)
Biocontainer bag distribution tower weight (empty)	440 kg (970 lb)
Bottle distribution tower dimensions (W × D × H)	210 × 70 × 133 cm (83 × 27 × 52 in.)
Bottle distribution tower weight (empty)	350 kg (772 lb)
Main system floor clearance	115 mm (4.5 in.)
Environmental conditions	5 – 30 °C, relative humidity 10 – 70% (non-condensing)
Materials of construction	Stainless steel 304
Surface finish	Cold rolled steel (typically Ra < 1 µm)
Ingress protection rating	IP54 (main, outer panel or chassis)

<sup>2</sup> Main system dimensions exclude the weighing platform. Weighing platform forms part of workstation dimensions.

## Utility specifications

Utilities	Specification
Electrical supply	230 VAC – 50 Hz (European), 208 VAC – 60 Hz (North American)
Operating voltage control	24 VDC – 20 A
Ingress protection for internal electrical cabinets	IP 56 (European), NEMA 4X (North American)
Recommended FI switch	RCD 30 mA, Type B
Power consumption	≤ 2.4 kVA
Amperage	10 A
Motor power	300 W (Primary pump), 140 W (Secondary pump)
Process air	Minimum pressure = highest integrity/leak test pressure + 1 bar , Maximum pressure = 10 bar, manual pressure regulator, pressure safety valve and Palltronic Flowstar IV Integrity Test Instrument
Instrument air	6 bar minimum, 10 bar maximum, 40 m <sup>3</sup> /h minimum flow rate, instrument air, clean, dry and oil free, pressure reducer and ON/OFF valve

## Component specifications

Component specifications, and sensor range and accuracies are as per OEM datasheets and correct at the time of compiling this proposal. Pall Corporation does not accept any responsibility in the case of deviation from the specifications outlined below.

Process equipment	Type	Specification
Primary pump	Watson-Marlow <sup>•</sup> 620R pump head	Flow: 0.15 – 10 L/min
Secondary pump	Watson Marlow Drivesure <sup>•</sup> 520R2	Flow: 0.12 – 3.5 L/min
Process valves	Gemü <sup>•</sup> Q30	Pneumatically operated pinch valve
Pressure	PendoTECH <sup>•</sup> – single-use sensor with pressure sensitivity chip	Range: -0.48 to 5.2 bar Accuracy: ± 2% from 0 to 0.41 bar, ± 3% from 0.41 to 2.07 bar, ± 5% from 2.07 to 4.1 bar
Level detector	Rechner <sup>•</sup> – capacitive sensor	Operating distance: 0.5 mm (min.), – 15 mm (max.)
Recovery tote	Plastic tray to support 5 L 2D biocontainer	High density polyethylene (HDPE)
Printer	Zebra <sup>•</sup> ZD420c – HC	Thermal transfer, 203 dpi, 15 – 108 mm label width
Handheld scanner	Zebra DS3678-HD3RF50BS1W	Wireless, Bluetooth <sup>•</sup> Class 1, image sensor 1280 × 800 pixels
Weighing platform	Mettler Toledo <sup>•</sup> PFK989 C600 Floor Platform	600 kg (max. capacity), recommended readability 1 g Accuracy: ± 0.5% at 1 kg <sup>1</sup> , ± 0.001% at 600 kg
Integrity test instrument	Palltronic Flowstar IV Integrity Test Instrument	Forward flow test: 0.1 – 1000 mL/ min., ± 3% of value or ± 0.05 mL/min, whichever is the greater

## Recommended process equipment and fittings

Process equipment	Recommendation
Product inlet mixer	LevMixer Stainless Steel Cubical Tank (ASME available upon request)
Product inlet mixer drive	LevMixer Drive Unit IV (LMG403)
Buffer flush feed tote	Allegro 100 L HDPE Tote
Buffer flush waste tote	Allegro 50 L Plastic 3D Tote
Bioprocessing workstation	Allegro Bioprocessing Workstation (LGRTTBSC)
Intermediate mixer	Allegro 50 L Standard Mixer and Weigh Platform

<sup>1</sup> Please note that the container workstation is ≥ 100 kg

## Automation specification

The system is equipped with Wonderware® Batch Management and Wonderware System Platform, which enables you to create recipes based on your individual process needs. Recipes are created using operations that are comprised of predefined phases. Each phase is configured to allow you to tailor the recipes as per your requirements.

The system also provides flexible phases, that allow for a fully manual creation of recipes.

Automation architecture is based on either Siemens S7 PLC or Rockwell CompactLogix PLC, an industrial PC, and a 22 in. touch screen HMI.

The automation platform allows the end user to create an unlimited number of recipes.

An OPC UA client is installed if a communication link is desired with the unit operation. The Pall Corporation automation platform enables compliance with 21 CFR Part 11 and follows the GAMP V life cycle for software development.

## Automation options

- PLC (Rockwell or Siemens) and HMI for local stand-alone control
- Remote I/O (no PLC) for integration into a DCS or SCADA system
- Remote I/O (no PLC) controlled by centralized PLC system


## Automation hardware specification<sup>1</sup>

Utilities	Specification
PLC vendor (option 1)	Siemens
PLC model (option 1)	SIMATIC® S7-1500, CPU 1516-3 PN/DP
PLC software (option 1)	Siemens TIA V16
PLC vendor (option 2)	Rockwell
PLC model (option 2)	5069-L380ERM
PLC software (option 2)	Rockwell Studio 5000 V32
Database	SQL
Operating system	Windows® Server 2016
Batch engine	Wonderware Batch Management
Historian	Wonderware Historian
Reporting	Dream Reports / Wonderware System Platform
HMI	SIMATIC ITC2200 V3 PRO, Industrial Thin Client, 22
Panel PC	INDUSTRY PC Stratus ztC Edge 200i
UPS	QUINT-UPS®/ 24DC
Remote I/O	Turck BL20® series
Ethernet switches	Phoenix Contact® 8 port (managed switch)
OPC server	Archestra® Client

<sup>1</sup> Technical equivalent replacements are possible depending on market availability, refer to Pall's Turnover Package (TOP) documentation.

## Process sequence

The chart below shows the separate process operations and process actions of the standard bulk fill process. Configurable phases are provided to cover these process steps. Examples are shown for both buffer and product-wet filter testing scenarios:

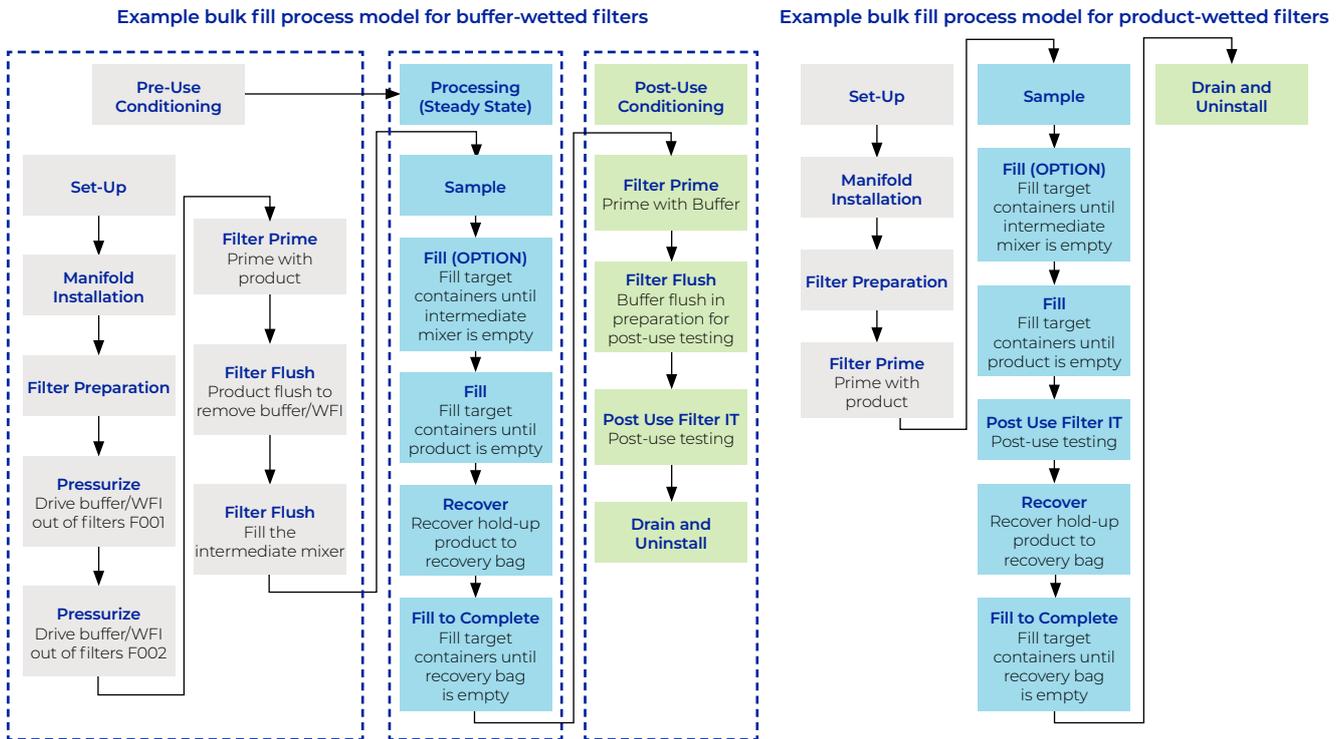

Additional phases which may be used to build custom recipes:




## Instructions for installation screens

The IFI have been created and are accessible via the HMI screen, providing you with a step-by-step guide to installing the single-use flow paths and making the relevant fluid connections. Sample IFI screens can be seen below:



## Process screens

Process screens have been created to summarize and expand on critical process information throughout the operation. The Palltronic Flowstar IV Integrity Test Instrument will also be controlled via the system HMI when performing filter integrity testing and/or pre-use manifold leak testing – sample screens of this procedure are also shown below.

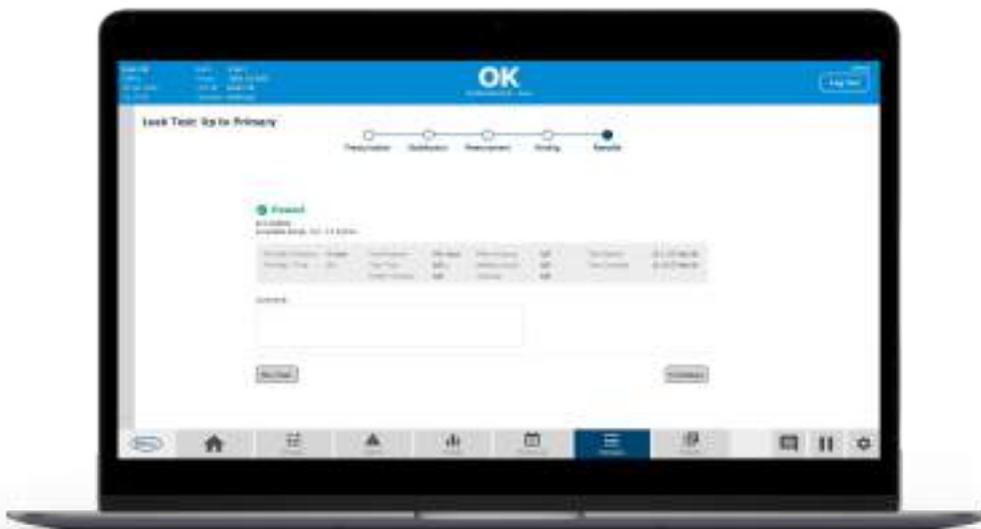

## Batch report

Upon operator selection, batch reports will be generated automatically at the end of a batch for each single step. Both a summary and detailed batch report is generated, and the content of these batch reports is predefined. Batch reports can be configured to specific needs by the end user via Wonderware Dream Reports. Sample batch reports can be provided upon request.

Predefined batch records contain the following major information:

- General batch information, phase information and transition conditions
- Global and recipe parameters including controller configuration
- Audit trails excerpt
- Alarms summary
- List of single-use manifold details that were registered during recipe execution
- Sample container list (including sample weight)
- Filled container list (including product weight)
- Trends for differential pressure across filters
- Trends for input weights
- Trends for filling weights

# Ordering information

## Automated system, workstations and towers

Part Number	Description
<b>Main system</b>	
ACBFSEUPLC	Allegro Connect Bulk Fill System: PLC 230 VAC, software automation
ACBFSWHPLC	Allegro Connect Bulk Fill System: PLC 208 VAC, software automation
ACBFSEUIO	Allegro Connect Bulk Fill System: I/O 230 VAC, DCS-ready no automation
ACBFSWHIO	Allegro Connect Bulk Fill System: I/O 208 VAC DCS-ready no automation
<b>All-in-one valve extension (or) tower</b>	
ACBFSBCREXT	Allegro Connect Bulk Fill System 6-valve biocontainer bag extension
ACBFSBTLEXT	Allegro Connect Bulk Fill System 6-valve bottle extension
ACBFSBCRTWRCE	Allegro Connect Bulk Fill System 16-valve biocontainer bag tower CE
ACBFSBCRTWRUL	Allegro Connect Bulk Fill System 16-valve biocontainer bag tower UL
ACBFSBTLWRCE	Allegro Connect Bulk Fill System 16-valve bottle tower CE
ACBFSBTLWRUL	Allegro Connect Bulk Fill System 16-valve bottle tower UL
<b>Biocontainer bag (or) bottle workstation</b>	
ACBFS12BCRWKSN	Allegro Connect Bulk Fill System 12-biocontainer bag workstation
ACBFS16BCRWKSN	Allegro Connect Bulk Fill System 16-biocontainer bag workstation
ACBFSBTLWKSN	Allegro Connect Bulk Fill System bottle workstation
<b>Bottle trays/inlays</b>	
ACBFS1LSQTRAY6	Allegro Connect Bulk Fill System 6 × 1 L square bottle tray
ACBFS2LSQTRAY6	Allegro Connect Bulk Fill System 6 × 2 L square bottle tray
ACBFS5LSQTRAY6	Allegro Connect Bulk Fill System 6 × 5 L square bottle tray
ACBFS10LSQTRAY6	Allegro Connect Bulk Fill System 6 × 10 L square bottle tray
ACBFS20LSQTRAY6	Allegro Connect Bulk Fill System 6 × 20 L square bottle tray
ACBFS1LSQTRAY16	Allegro Connect Bulk Fill System 16 × 1 L square bottle tray
ACBFS2LSQTRAY16	Allegro Connect Bulk Fill System 16 × 2 L square bottle tray
ACBFS5LSQTRAY16	Allegro Connect Bulk Fill System 16 × 5 L square bottle tray
ACBFS10LSQTRAY12	Allegro Connect Bulk Fill System 12 × 10 L square bottle tray
ACBFS20LSQTRAY12	Allegro Connect Bulk Fill System 12 × 20 L square bottle tray
<b>Seismic feet and UL certification</b>	
ACSEISMIC	Allegro Connect system seismic leveling feet
ACSEISMICBCRTWR	Allegro Connect system seismic leveling biocontainer bag tower
ACSEISMICBTLWR	Allegro Connect system seismic leveling bottle tower
ACULCERT	Allegro Connect equipment UL certificate 61010
<b>FAT</b>	
ACBFSFAT	Allegro Connect Bulk Fill System: FAT 2.5 days
ACBFSFATEXT	Allegro Connect Bulk Fill System: Extended FAT 5 days
<b>Other accessories</b>	
ACGEN3COMSCBL	Allegro Connect system GEN3 mixer communication cable
ACGEN4COMSCBL	Allegro Connect system GEN4 mixer communication cable
ACLGRCOMSCBL	Allegro Connect system LGR mixer communication cable

## Ordering guide suggestion for system hardware



All-in-one biocontainer bags



All-in-one bottles



Tower biocontainer bags



Tower bottles



All-in-one hybrid\*  
(biocontainer bags and bottles)

\* Towers can also be used with the hybrid system

### Step 1: Choose the central unit (with or without automation)<sup>1</sup>

Central bulk fill unit for CE region with automation	Central bulk fill unit for CE region without automation	Central bulk fill unit for UL region with automation	Central bulk fill unit for UL region without automation
ACBFSCEPLC	ACBFSCEIO	ACBFSULPLC	ACBFSULIO

<sup>1</sup> 1 × label roll and 1 × printer cartridge are included with the system. For additional quantity please place the order with Codeway Ltd (England) for vendor P/N: 8000T CryoCool (labels) and 4800 Resin (cartridge)

### Step 2: Choose the system configuration for biocontainer bags/bottles/hybrid based on container size and number of fills

All-in-one biocontainer bag(s) for 1 L to 20 L container size up to 120 fills	All-in-one bottle for 1 L to 20 L container size up to 120 fills	All-in-one hybrid (biocontainer bag(s) and bottle) for 1 L to 20 L container size up to 120 fills	Bottle distribution tower for 1 L to 5 L container size from 121 to 320 fills
ACBFSBCREXT	ACBFSBTLEXT	ACBFSBCREXT and ACBFSBTLEXT	ACBFSBTLTWRCE/ ACBFSBTLTWRUL
ACBFSPFK9C600	ACBFSPFK9C600	ACBFSPFK9C600	ACBFSPFK9C600
ACBFS12BCRWKSN	ACBFSBTLWKSN <sup>2</sup>	ACBFS12BCRWKSN and ACBFSBTLWKSN <sup>2</sup>	ACBFSBTLWKSN <sup>2</sup>
Bottle distribution tower for 10 L and 20 L container size from 121 to 240 fills	Biocontainer bag(s) distribution tower for 1 L to 5 L container size from 121 to 320 fills	Biocontainer bag(s) distribution tower for 10 L and 20 L container size from 121 to 320 fills	
ACBFSBTLTWRCE/ ACBFSBTLTWRUL	ACBFSBCRTWRCE/ ACBFSBCRTWRUL	ACBFSBCRTWRCE/ ACBFSBCRTWRUL	
2 × ACBFSPFK9C600	ACBFSPFK9C600	2 × ACBFSPFK9C600	
2 × ACBFSBTLWKSN <sup>2</sup>	ACBFS16BCRWKSN <sup>2</sup>	2 × ACBFS16BCRWKSN <sup>2</sup>	

<sup>2</sup> Proceed to step 2A to choose complementary trays/inlays. Step 2A and Step 3 are only applicable for bottle filling configuration.



Step 2A: Choose 2 complementary single bottle trays (inlays) or 1 complementary dual bottle tray/inlay for each ACBFSBTLWKS depending on the size of the bottles used in the process

Available sizes	Part number	Available sizes	Part number
<b>Single bottle tray/inlay</b>		<b>Dual bottle tray/inlay</b>	
6 × 1 L	ACBFS1LSQTRAY6	16 × 5 L	ACBFS5LSQTRAY16
6 × 2 L	ACBFS2LSQTRAY6	12 × 10 L	ACBFS10LSQTRAY12
6 × 5 L	ACBFS5LSQTRAY6	12 × 20 L	ACBFS20LSQTRAY12
6 × 10 L	ACBFS10LSQTRAY6	N/A	N/A
6 × 20 L	ACBFS20LSQTRAY6	N/A	N/A
16 × 1 L	ACBFS1LSQTRAY16	N/A	N/A
16 × 2 L	ACBFS2LSQTRAY16	N/A	N/A

Step 3 (optional): Choose additional bottle trays/inlays as required

Available sizes	Part number	Available sizes	Part number
<b>Single bottle tray/inlay</b>		<b>Dual bottle tray/inlay</b>	
6 × 1 L	ACBFS1LSQTRAY6	16 × 5 L	ACBFS5LSQTRAY16
6 × 2 L	ACBFS2LSQTRAY6	12 × 10 L	ACBFS10LSQTRAY12
6 × 5 L	ACBFS5LSQTRAY6	12 × 20 L	ACBFS20LSQTRAY12
6 × 10 L	ACBFS10LSQTRAY6	N/A	N/A
6 × 20 L	ACBFS20LSQTRAY6	N/A	N/A
16 × 1 L	ACBFS1LSQTRAY16	N/A	N/A
16 × 2 L	ACBFS2LSQTRAY16	N/A	N/A

### Single-use assemblies

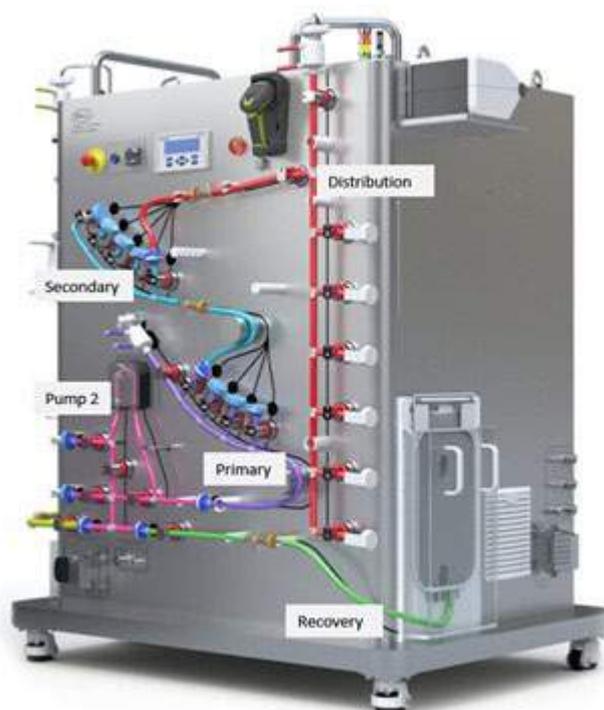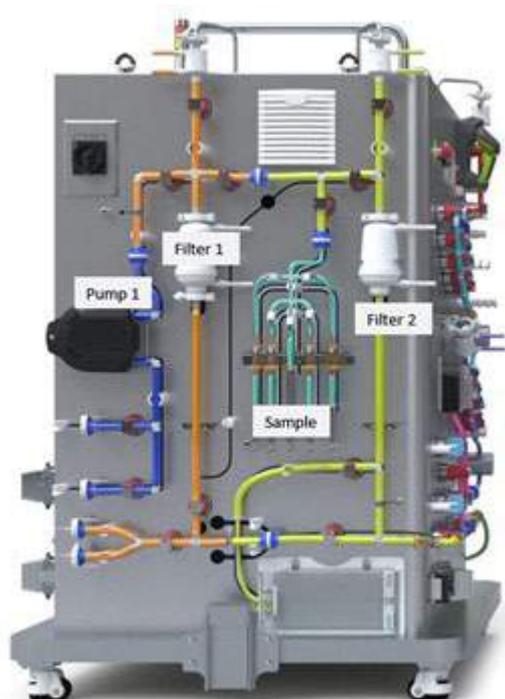

## Part number matrix

The flow path selection table shown below lists the standard part numbers of the Pall products designed and approved to support the Allegro Connect Bulk Fill System. Due to the configurability of the system enabling numerous combinations of the filter capsules (membrane and size), we are only able to list a limited number of the possible filtration assemblies in the table below. Filtration combination sets, which do not appear in this table but are within the scope of the system, are available from Pall Corporation, on request.

Flow path selection	Part number	Description
<b>Step 1: choose pump 1 manifold</b>		
Pump 1	9496-1356Z	Pump 1 manifold with Kleenpak Presto Sterile Connector
<b>Step 2: choose what type of filtration manifold is required for the process</b>		
<b>No filtration</b>		
Bypass manifold	9496-1579R	Filter interlink manifold with Kleenpak Presto Sterile Connector
Sampling manifold	9496-1376Z	Sampling manifold with 4 x 50 mL biocontainer bags
	9496-1579S	Serial 2 manifold with no filter
<b>Single filtration (select bypass and sampling manifold. Select a serial 2 filter)</b>		
Bypass manifold	9496-1579R	Filter interlink manifold with Kleenpak Presto Sterile Connector
Sampling manifold	9496-1376Z	Sampling manifold with 4 x 50 mL biocontainer bags
	9496-1572T	Serial 2 manifold with KA1EKVP16G filter
	9496-1553W	Serial 2 manifold with NP1LUECVPIG filter
	9496-1561U	Serial 2 manifold with NP5LUECVPI6G filter
	9496-1579K	Serial 2 manifold with NP6UECVPIG filter
	9496-1358F	Serial 2 manifold with NP7EKVP16G filter
	9496-1581F	Serial 2 manifold with NP7UECVPIG filter
	9496-1800S	Serial 2 manifold with KA02EKVP1G filter
	9496-1785E	Serial 2 manifold with KA3EKVP16G filter
<b>Serial filtration (select sampling manifold. Select a serial 1 and serial 2 combination)</b>		
Sampling manifold	9496-1376Z	Sampling manifold with 4 x 50 mL biocontainer bags
	9496-1558Q	Serial 1 manifold with KA1EKVP16G filter
	9496-1572T	Serial 2 manifold with KA1EKVP16G filter
	9496-1519F	Serial 1 manifold with KA3EAVPIG filter
	9496-1553W	Serial 2 manifold with NP1LUECVPIG filter
	9496-1557E	Serial 1 manifold with NP5LUEAVPIG filter
	9496-1561U	Serial 2 manifold with NP5LUECVPI6G filter
	9496-1579J	Serial 1 manifold with NP6UEAVP6G filter
	9496-1579K	Serial 2 manifold with NP6UECVPIG filter
	9496-1357E	Serial 1 manifold with NP7UEAVPIG filter
	9496-1358F	Serial 2 manifold with NP7EKVP16G filter
<b>Parallel filtration (select sampling manifold. Select a parallel 1 and parallel 2 combination)</b>		
Sampling manifold	9496-1376Z	Sampling manifold with 4 x 50 mL bags
	9496-1574M	Parallel 1 manifold with KA1EKVP16G filter
	9496-1574P	Parallel 2 manifold with KA1EKVP16G filter
	9496-1560Z	Parallel 1 manifold with NP5LUECVPI6G filter
	9496-1560Y	Parallel 2 manifold with NP5LUECVPI6G filter
	9496-1579M	Parallel 1 manifold with NP6EKVP6G filter

	9496-1579N	Parallel 2 manifold with NP6EKVP6G filter
	9496-1581H	Parallel 1 manifold with NP7UECVPIG filter
	9496-1581G	Parallel 2 manifold with NP7UECVPIG filter
	9496-1358G	Parallel 1 manifold with NP7EKVP16G filter
	9496-1513S	Parallel 2 manifold with NP7EKVP16G filter
<b>Step 3: choose pump 2, recovery and multiplier manifolds (*quantity of secondary multiplier manifold varies based on number of containers to be filled)</b>		
Pump 2	9496-1362T	Pump 2 intermediate mixer manifold
Primary	9496-1577Z	Primary multiplier manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
Secondary (if required)*	9496-1578B	Secondary multiplier manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
Recovery bag	9496-1402Z	Recovery bag manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
<b>Step 4: choose distribution manifolds (quantity of distribution manifold varies based on number of containers to be filled)</b>		
Biocontainer bag distribution ('all-in-one')	9496-1579F	6 x biocontainer bag distribution manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
Bottle distribution ('all-in-one')	9496-1519E	6 x bottle distribution manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
Biocontainer bag tower distribution ( $\leq$ 5L)	9496-1579G	16 x biocontainer bag distribution manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
Biocontainer bag tower distribution ( $>$ 5L)	9496-1579T	12 x biocontainer bag distribution manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
Bottle tower distribution ( $\leq$ 5L) <sup>1</sup>	9496-1513T	16 x bottle distribution manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
Bottle tower distribution ( $>$ 5L)	9496-1519D	12 x bottle distribution manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
<b>Step 5: choose storage biocontainer bag(s) or bottle cap manifold (quantity of storage containers varies based on the number of containers to be filled)</b>		
<b>Biocontainer bag(s)</b>		
	7190-1561S	1 L storage bag manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
	7190-1561W	2 L storage bag manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
	7190-1561X	5 L storage bag manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
	7190-1561T	10 L storage bag manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
	7190-1571R	20 L storage bag manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter
<b>Bottle cap</b>		
Bottle cap manifold for tower	7496-1579U	Bottle cap manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter (for tower)
Bottle cap manifold for all-in-one	7496-1671Y	All-in-one bottle cap manifold with Kleenpak Presto Sterile Connector and Kleenpak Sterile Disconnecter

<sup>1</sup> Manifold can be used for 16 x filling with 10 L and 20 L bags when additional weighing scale and 16-biocontainer bag workstation is used

In order to reduce our carbon footprint, we strive to provide single-use systems manufactured regionally. However, to ensure security of supply you may receive product from multiple global sites.

## Scientific and laboratory services

The scientific and regulatory knowledge that supports the selection, adoption, and ongoing use of critical process technology – coupled with analytical, imaging and measurement capabilities – creates a versatile and practical resource ready to respond to an ever-changing industry. Pall duplicates these laboratories across the globe and leverages their cumulative knowledge to deliver practical scientific and regulatory support to all process technologies, to keep you moving forward.

## Technical services

The accessibility of local technical support networks minimizes delays in your journey at all points. From the early stage of process development to on-site support for mature processes, Pall's technical support groups are there to help remove barriers to progress and make your journey as rapid and stress-free as possible. Our knowledge of the technology and the process can be applied to everything from training to troubleshooting and consultancy. Our global team of technology experts is on hand to respond to your changing needs.

## Advanced separation systems

Operating within the defined design space demands the monitoring and control of critical process parameters to assure product quality. Systems that control critical unit operations and communicate with your existing process components can control process risks and maximize productivity by reducing operator involvement for many processes. Pall applies strong engineering and regulatory understanding to deliver compliant and qualified systems that safeguard and simplify your journey.

## Process development services

Prior knowledge is a rare and valuable commodity, especially when preparing to take a new direction or when under pressure to deliver to a tight deadline. Take advantage of Pall's experience, process knowledge and technical know-how to help you achieve your goals. From the optimization of an end-to-end continuous process to establishing the right parameters for a single unit operation, our teams of scientists are ready to work with you and to generate the data you need to make the critical decisions necessary for success.

## Validation services

Arriving at your destination counts for nothing without the necessary paperwork to proceed to the next stage. Pall's Validation Services are committed to delivering the supporting data packages and analysis required to quantify process risk and support regulatory submission. Our strengths include critical filtration technologies, such as the performance validation of sterilizing grade filtration, and we are at the forefront of the evolving needs in the area of extractables and leachables for all product contact components. We combine the generation of data with interpretation and consultancy to deliver data packages that are ready for regulatory scrutiny, and to ensure there are no barriers to progress.

## Servicing and maintenance

Our range of service packages keeps your equipment protected and well maintained, and includes itemized, pay-as-you-go services, start-up care and training packages, and a variety of post-warranty service plans that include priority response times, discounts for emergency repairs, and flexible payment options. Pall service plans provide total peace of mind and worry-free support throughout the coverage period.



Corporate Headquarters  
Port Washington, NY, USA  
+1-800-717-7255 toll free (USA)  
+1-516-484-5400 phone

European Headquarters  
Fribourg, Switzerland  
+41 (0)26 350 53 00 phone

Asia-Pacific Headquarters  
Singapore  
+65 6389 6500 phone

Visit us on the Web at [www.pall.com/biotech](http://www.pall.com/biotech)  
Contact us at [www.pall.com/contact](http://www.pall.com/contact)

Pall Corporation has offices and plants throughout the world. To locate the Pall office or distributor nearest you, visit [www.pall.com/contact](http://www.pall.com/contact).

The information provided in this literature was reviewed for accuracy at the time of publication. Product data may be subject to change without notice. For current information consult your local Pall distributor or contact Pall directly.

© Copyright 2022, Pall Corporation. Pall, , Allegro, Kleenpak, LevMixer and Palltronic are trademarks of Pall Corporation. ® Indicates a trademark registered in the USA.

◆ ArchestrA is a trademark of Aveva Software LLC, Aspen is a trademark of The Aspen Institute, Bluetooth is a registered trademark of The Bluetooth Special Interest Group (SIG), CompactLogix is a trademark of Rockwell International Corporation, CryoCool is a trademark of Thermo Fisher Scientific Inc., Drivesure is a trademark of Watson-Marlow Ltd, GAMP is a trademark of the International Society for Pharmaceutical Engineering, Gemü is a trademark of the GEMÜ Group, Mettler Toledo is a trademark of Mettler-Toledo GmbH, PendoTECH is a trademark of PendoTECH, Phoenix Contact is a trademark of Phoenix Contact GmbH & Co. KG, QUINT-UPS is a trademark of Quint Holding B.V., Rechner is a trademark of Rechner Industrie-Elektronik GmbH, Rockwell is a trademark of Rockwell International Corporation, Siemens is a trademark of Siemens AG, SIMATIC is a trademark of Siemens AG, Turck BL20 is a trademark of Hans Turck GmbH & Co. KG, Watson-Marlow is a trademark of Watson-Marlow Ltd, Windows is a trademark of Microsoft Corporation, Wonderware is a trademark of Aveva Software LLC.

USD3665/22.7985  
06/2022